


Protocol for IHC-Fr(Anti Iba, Rabbit)

Wistar Rat (7 week old, ♂) was perfusion-fixed with 4% paraformaldehyde. replaced sucrose, and prepared 50µm brain section by microtome.

- Wash :** 0.3 % TritonX-100 in PBS、 5 min×3
↓
Blocking : 1 % BSA, 0.3 % TritonX-100/PBS, 2hour , RT
↓
Primary antibody : rabbit anti-Iba1(1 : 1000), 1 % BSA, and 0.3 % TritonX-100 in PBS、
Over night 、 4°C
↓
Wash : 0.3 % TritonX-100 in PBS、 5 min×3
↓
Secondary antibody : AlexaFluor488 anti-rabbit IgG(1:1000), 1 % BSA, and 0.3 %
TritonX-100 in PBS, 2hour, RT
↓
Wash : 0.3 % TritonX-100 in PBS、 5 min×3
↓
Mount

IHC data sample

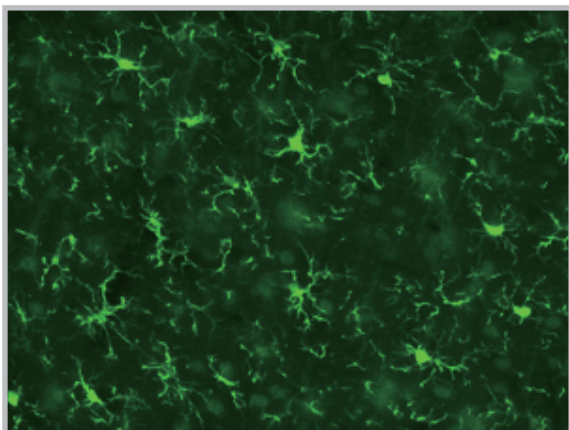

产品列表

产品编号	产品名称	包装
019-19741	小胶质细胞/巨噬细胞特异性蛋白抗体 (免疫组化) Anti Iba1, Rabbit (for Immunocytochemistry)	50 µg

上述试剂仅供实验研究用,不可用作“医药品”、“食品”、“临床诊断”等。

Listed products are intended for laboratory research use only, and not to be used for drug, food or human use. / Please visit our online catalog to search for other products from FUJIFILM Wako; <https://labchem-wako.fujifilm.com> / This leaflet may contain products that cannot be exported to your country due to regulations. / Bulk quote requests for some products are welcomed. Please contact us.

富士胶片和光(广州)贸易有限公司

www.boppard.cn

广州市越秀区先烈中路69号东山广场30楼
3002-3003室
询价: info@boppard.cn
目录价查询: bb-china.net

北京 Tel: 010 64136388
上海 Tel: 021 62884751
广州 Tel: 020 87326381
香港 Tel: 852 27999019

官方微信



目录价查询

