

# IHC Protocol (Paraffin Section) Anti Iba1, Rabbit

100% xylene, 5 min × 2  
100% Ethanol, 5 min × 2  
0.3% H<sub>2</sub>O<sub>2</sub> in Methanol, 20 min, (inactivation of endogenous peroxidase)  
90% Ethanol, 2 min  
70% Ethanol, 2 min  
H<sub>2</sub>O 3 min  
PBS 5 min × 2  
↓  
Citrate Buffer (pH6.0) 90°C, 9 min  
Incubate to RT  
↓  
PBS 5 min × 2  
↓  
2% BSA in PBS RT, 60min (blocking)  
↓  
PBS 5 min × 2  
↓  
Anti Iba1, Rabbit (1:1,000) O/N  
↓  
PBS 5 min × 2  
↓  
Secondary antibody (Biotinylated anti Rabbit IgG), 60 min, RT  
↓  
PBS 5 min × 2  
↓  
ABC, 1 hour, RT  
↓  
PBS, 5 min × 2  
↓  
DAB solution  
↓  
PBS, 5 min × 2  
↓  
Mount

## 产品列表

产品编号	产品名称	包装
019-19741	小胶质细胞/巨噬细胞特异性蛋白抗体(免疫组化) Anti Iba1, Rabbit (for Immunocytochemistry)	50 µg

上述试剂仅供实验研究用,不可用作“医药品”、“食品”、“临床诊断”等。

Listed products are intended for laboratory research use only, and not to be used for drug, food or human use. / Please visit our online catalog to search for other products from FUJIFILM Wako; <https://labchem-wako.fujifilm.com> / This leaflet may contain products that cannot be exported to your country due to regulations. / Bulk quote requests for some products are welcomed. Please contact us.

富士胶片和光(广州)贸易有限公司

[www.boppard.cn](http://www.boppard.cn)

广州市越秀区先烈中路69号东山广场30楼  
3002-3003室  
询价: [info@boppard.cn](mailto:info@boppard.cn)  
目录价查询: [bb-china.net](http://bb-china.net)

北京 Tel: 010 64136388  
上海 Tel: 021 62884751  
广州 Tel: 020 87326381  
香港 Tel: 852 27999019

官方微信



目录价查询

